

Standard operating procedure of Micro centrifuge

- 1.Objective: This instruction explains how use and keep the micro centrifuge apparatus
- 2.Scope: using for centrifuging the different ranges of solutions with maximum 2 mL volume.
- 3.Responsibility: Laboratory assistance is responsible for the accuracy of the apparatus and should be reported to him/her in case of problems
- 4.Materials and equipment:...
- 5.Guidelines for use :
 - Turn on the apparatus
 - Open the door of apparatus with blue key
 - Place the tubes in face to face pattern for maintaining the balance of apparatus
 - Adjustment the time and centrifuging speed of the apparatus with the blue set bottoms and run your reaction.
- 6.Warning and safety precautions:
If any abnormal sound or vibrate are felt, press the “Stop” bottom immediately .
- 7.Maintenance :
The door of the apparatus should be always closed.

Document ID: **CoreLab. 15**

Document name	Standard operating procedure of Micro centrifuge
Date	99/4/1
Full file name	Standard operating procedure of Micro centrifuge
Document description	These instructions describe the usage, maintenance and calibration of Micro centrifuge which is kept in Core laboratory.
Producer	Core Lab
